

**Formula \* in g/L**

|                                     |                       |
|-------------------------------------|-----------------------|
| Enzymatic digest of casein.....     | 10.000                |
| Yeast extract.....                  | 1.000                 |
| Magnesium chloride (anhydrous)..... | 28.100 <sup>(*)</sup> |
| Sodium chloride .....               | 5.000                 |
| Malachite green.....                | 0.010                 |
| Potassium Chlorate.....             | 1.000                 |

Final pH 6.9 ±0,2 at 25 °C

(\*) Equivalent to 60.0 g of MgCl<sub>2</sub>·6H<sub>2</sub>O

\* Adjusted and /or supplemented as required to meet performance criteria

**Directions**

Dissolve 45.11 g of the powder in 1 liter of distilled water and sterilize in the autoclave for 15 minutes at 121 ° C. It is allowed to cool to 50 °C and add 1 ml of a sterile solution of Ticarcillin and Triclosan (Irgasan®) at 0,1%. Homogenize and distribute in suitable containers trying to minimize contact with air (relative anaerobiosis).

**Description**

This medium was originally formulated by Wauter *et al.*, starting from the enrichment broth for salmonellae from Rappaport *et al.*, to which they modified the proportions of magnesium chloride and malachite green and added potassium chlorate that inhibits the growth of nitratase A-producing enterobacteria. The selectivity for *Yersinia enterocolitica* is achieved with the inclusion of triclosan (Irgasan®) that acts against gram-positive microorganisms and ticarcillin, an antibiotic that interferes with the formation of bacterial cell walls.

This medium works very well to detect the *Y. enterocolitica* pathogens of biotype 4 and serotype O:3, but it is not suitable for detecting other serotypes.

**Technique**

Proceed according to current national or international standards, established and tested protocols or according to the procedures established and accepted in each laboratory.

**Quality control**

**Incubation temperature:** 25±1 °C

**Incubation time:** 44± 4h

**Inoculum:** Practical range 50 - 100 CFU (Productivity)/ 10<sup>4</sup>-10<sup>6</sup> CFU (Selectivity) according to ISO 11133:2014/ Amd 1:2018

**Microorganism**
**Growth**
**Remarks**

|  |                          |                      |
|--|--------------------------|----------------------|
| <i>Yersinia enterocolitica</i> DSM® 13030 +(1)+(2) | Good                     | > 10 CFU in CIN Agar |
| <i>Escherichia coli</i> ATCC® 25922 (1)            | Total Inhibited          | -                    |
| <i>Pseudomonas aeruginosa</i> ATCC® 27853 (2)      | Total Inhibited          | -                    |
| <i>Proteus mirabilis</i> ATCC® 29906               | Partial/ total Inhibited | <10 CFU in TSA       |

---

### References

- ISO Standard 10273 (2017) Microbiology of food and animal feeding stuffs - Horizontal method for the detection of presumptive pathogenic *Yersinia enterocolitica*.
- ACUFF, G.R. (2001) Media, Reagents and Stains in Compendium of methods of microbiological examination of foods. 3<sup>rd</sup> E d. Chap. 62. APHA. Washington D.C.
- ATLAS, R.M. & L.C. PARKS (1993) Handbook of Microbiological Media. CRC Press, Inc. London
- CORRY, J.E.L., G.D.W. CURTIS & R.M. BAIRD (2012) Handbook of culture media for food and water microbiology. RSC Publishing. Cambridge. UK.
- CURTIS, G.D.W. & R.M. BAIRD (1993) Pharmacopoeia of Culture Media for Food Microbiology: Additional monographs (II). Int. J. Food Microbiol 17:201-266
- de BOER, E. (2012) Culture media for the isolation of *Yersinia enterocolitica* from foods. Chap 15 in Handbook of culture media for food and water microbiology. (Corry *et al.*, eds) RSC Publishing. Cambridge. UK.
- de ZUTTER, L., L. LeMORT, M. JANSSENS & G. WAUTERS (1994) Short-comings of Irgasan Ticarcillin Chlorate Broth for the enrichment of *Yersinia enterocolitica* biotype 2 serotype 9 from meat. Int. J. Food Microbiol. 23:231-237
- de ZUTTER, L., M. JANSSENS & G. WAUTERS (1995) Detection of *Yersinia enterocolitica* serogroup O:3 using different inoculation methods of the enrichment médium Irgasan Ticarcillin Chlorate. Contrib. Microbiol. Immunol. 13:123-125
- ISO Standard 10273:2017 Microbiology of the food chain. Horizontal method for the detection of pathogenic *Yersinia enterocolitica*.
- DOWNES, F.P. & K. ITO (2006) Compendium of methods for the microbiological examination of foods. 4<sup>th</sup> Ed. APHA. Washington D.C.
- PRPIC, J.K & D. HUGHES (1989) *Yersinia enterocolitica* in Foodborne Microorganisms of Public Health Significance 4<sup>th</sup> Ed. (K.A. Buckle *et al.* Eds) Chap. 6. AIFST. Australia.
- RAPPAPORT, F., N. KONFORTIN & B. NAVON (1956) A new enrichment médium for certain *Salmonellae*. J. Clin. Pathol. 9:261-266
- SCHIEMANN, D.A. & G. WAUTERS (2001) *Yersinia* in Compendium of methods for the microbiological examination of foods. 3<sup>rd</sup> ed. Chap. 27. APHA. Washington D.C.
- VANDERZANT, C. & D. SPLITTERSTOESSER(Eds) (2001) Compendium of methods for the microbiological examination of foods. 3<sup>rd</sup> ed. APHA. Washington D.C.
- WAUTERS, G., V. GOOSENS, M. JANSSENS & J. VANDEPITTE (1988) New enrichment method for isolation of pathogenic *Yersinia enterocolitica* serogroup O:3 from pork. Appl. Environ. Microbiol. 54:851-854
- WEAGANT, S.D. & P. FENG (2006) *Yersinia* in Compendium of methods for the microbiological examination of foods. 4<sup>th</sup> ed. APHA. Washington D.C.

### Storage

For laboratory use only. Keep tightly closed, away from bright light, in a cool dry place (+4 °C to 30 °C).